

MICROLINKS TECHNOLOGY CO.,LTD

User's Manual

Microscope Installation Manual

MLTC

2014/05/22 Update Ver.2.8A

Content

- 1. **Please insert CD in CD-Rom**3
- 2. **Install and uninstall microscope software**.....3
 - 2.1 Install software.....4
 - 2.1.1 Setup software4
 - 2.1.2 Software Setup Wizard4
 - 2.1.3 Choose installation route5
 - 2.1.4 Installing.....5
 - 2.1.5 Install Completed6
 - 2.1.6 Software icon on desktop6
 - 2.2. Uninstall Microscope software7
 - 2.2.1 Choose to uninstall Microscope software.....7
 - 2.2.2 Confirm to Uninstall7
 - 2.2.3 Uninstalling.....7
 - 2.2.4 Uninstall Completed.....7

1. Please insert CD in CD-Rom

The package provides setup CD. If the setup CD is missing or damaged, please download setup files from <http://www.vitiny.com>. Please see the following steps to complete the Microscope installation.

- **Step 1: Please insert CD.**
- **Step 2: Install Microscope software. Please refer to section 2.1.**

To uninstall Microscope software in your PC, please follow below 2 steps to uninstall.

- **Step 1: Uninstall Microscope software. Please refer to section 2.2.**

The install window will pop out automatically once CD is inserted.



Fig 1-1 Install window

If the install window does not show, please click Autorun.exe from CD, the same install window will show up, too.



Fig 1-2 Autorun.exe

2. Install and uninstall microscope software

Please select the Fig.1-1 to start setup Microscope software program.

2.1 Install software

2.1.1 Setup software

First, click “Install Microscope software”, next, please select a language, see Fig. 2-1, the initial language will follow up the pc OS system. If you want to cancel setup please click “Cancel” then click “OK”

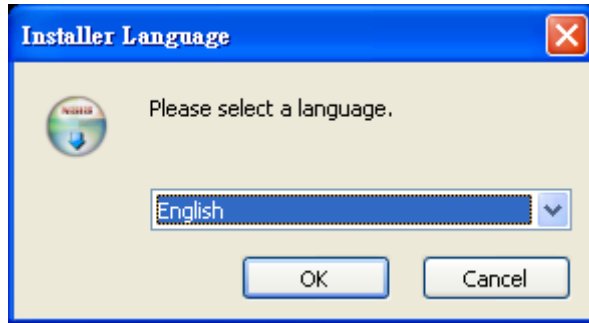


Fig 2-1seup language window

2.1.2 Software Setup Wizard

After choosing the language, go to welcome to the Microscope setup wizard window see Fig.2-2, then please click “Next” then continue. If want to cancel please click “Cancel” button. In the following figs, “name” represents Microscope’s software name, “X.XXX.XXXx” represents the software version.

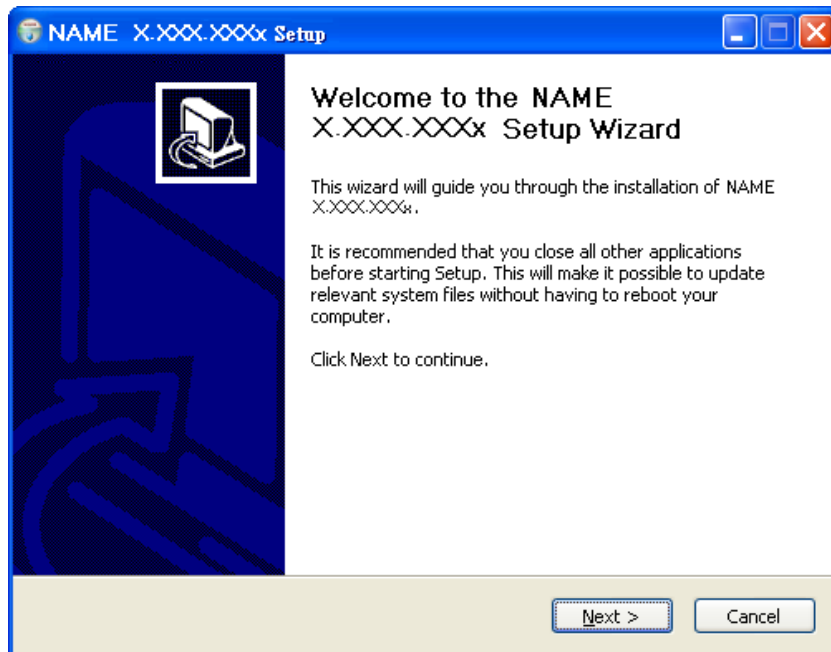


Fig 2-2 welcome to the Microscope setup wizard

2.1.3 Choose installation route

Next step please choose the installation route window see Fig.2-3. If want to change file route click “Browse” to change setup path, then please click “install” then start installing.

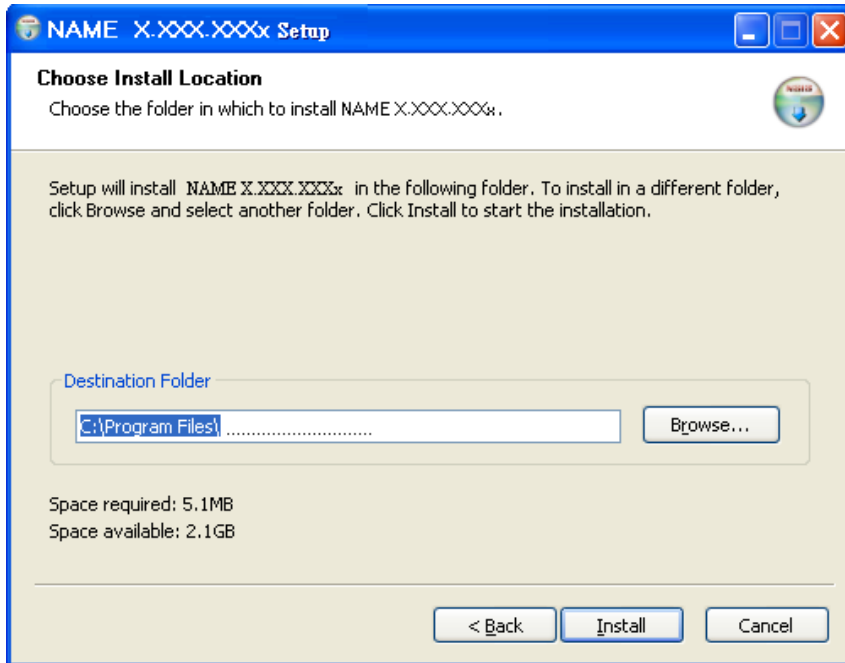


Fig 2-3 choose installation route

2.1.4 Installing

Continue install wizard will start installing program ad directory Fig 2-4.

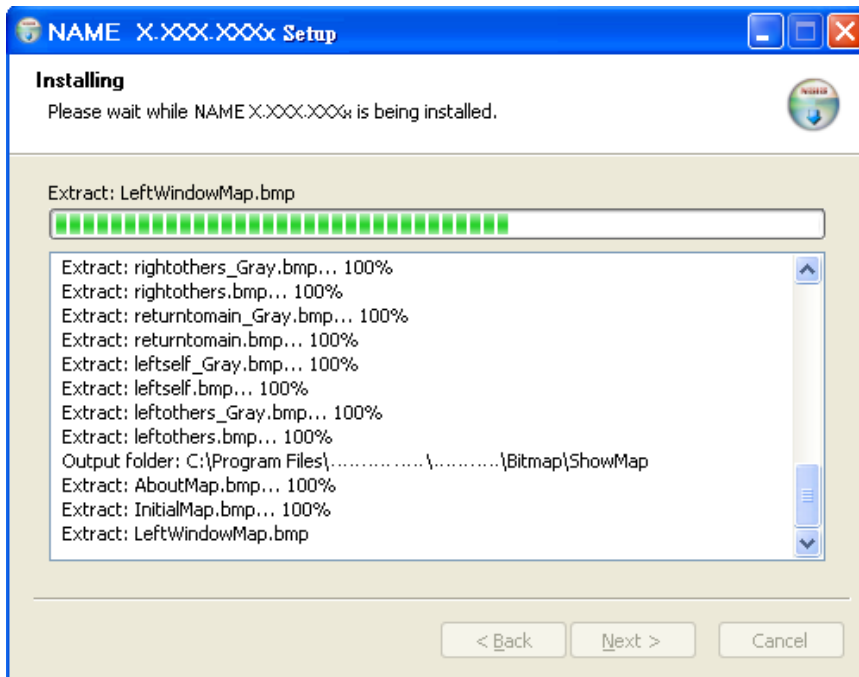


Fig 2-4 installing

2.1.5 Install Completed

When the Microscope program complete installation (Fig 2-5), please click “Finish”.

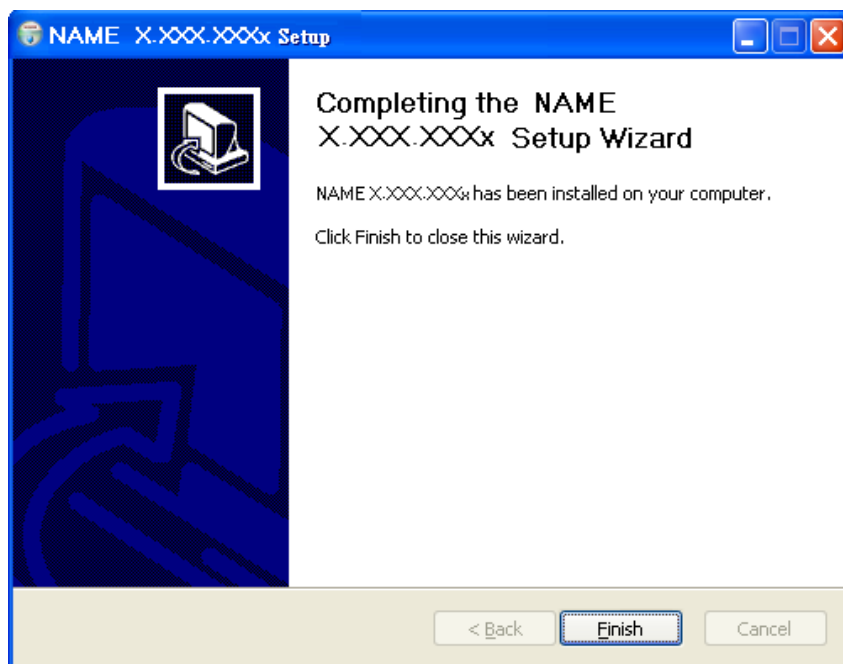


Fig 2-5 Microscope wizard install completed

2.1.6 Software icon on desktop

Microscope software shortcut will show on the Desktop as Fig.2-6.



Fig 2-6 Microscope shortcut icon

Microscope shortcut (Fig 2-6) will show on the Desktop. Or go “Start”-> All Programs-> MicroLinks Technology -> Microscope (see Fig 2-7).

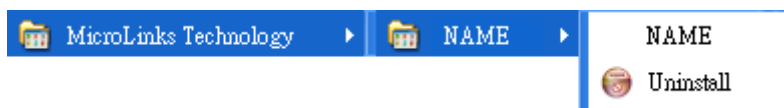


Fig 2-7 Microscope in All Programs

2.2. Uninstall Microscope software

2.2.1 Choose to uninstall Microscope software

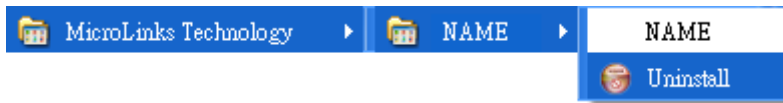


Fig.2-8 Microscope software Uninstall

2.2.2 Confirm to Uninstall

Please make sure if you want to completely remove Microscope software. See Fig.2-9

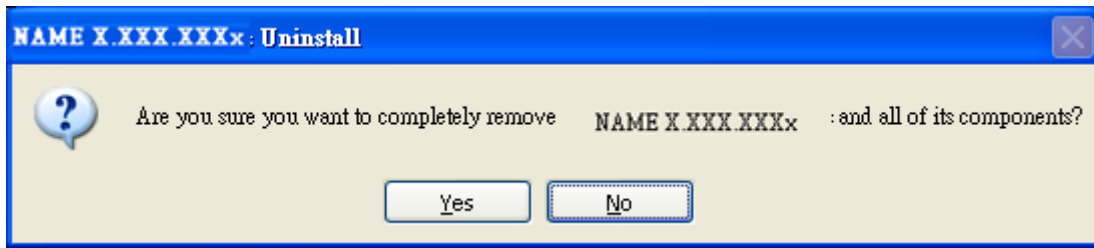


Fig.2-9

2.2.3 Uninstalling

In processing of uninstall the Microscope software, see Fig.2-10

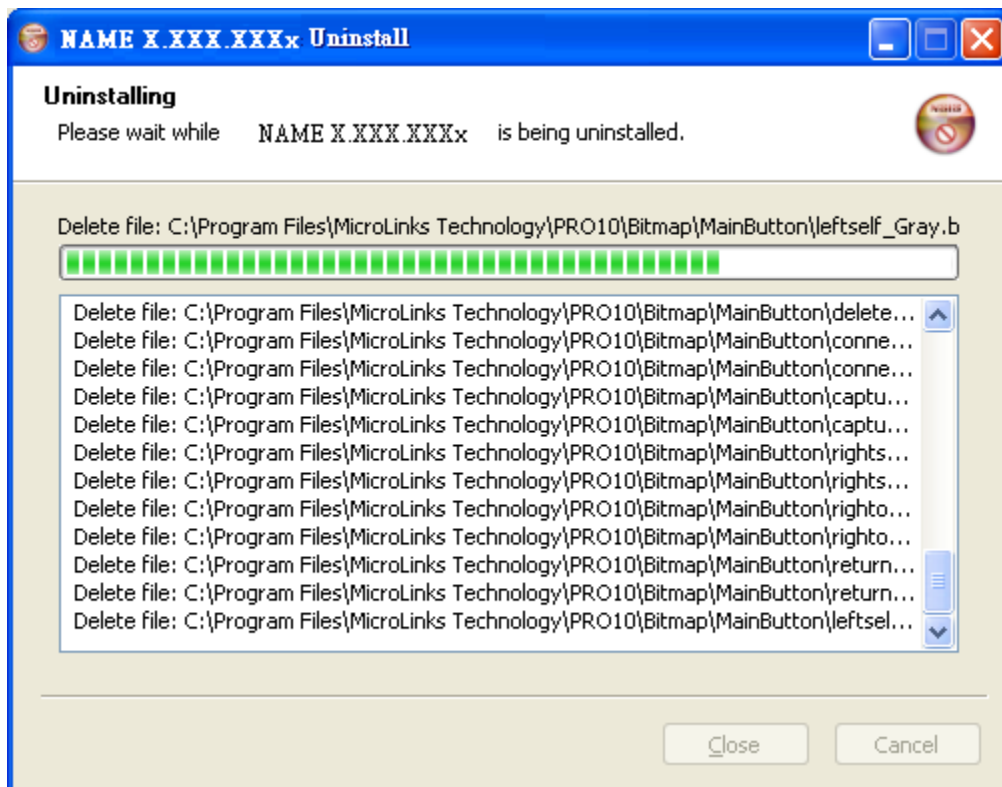


Fig.2-10 In process

2.2.4 Uninstall Completed

Microscope software was successfully removed from your PC, please click OK to finish.



Fig. 2-11 Uninstall Completed